



Additivity of Heats of Reaction: Hess's Law

In this experiment, you will use a Styrofoam-cup calorimeter to measure the heat released by three reactions. One of the reactions is the same as the combination of the other two reactions. Therefore, according to Hess's law, the heat of reaction of the one reaction should be equal to the sum of the heats of reaction for the other two. This concept is sometimes referred to as the *additivity of heats of reaction*. The primary objective of this experiment is to confirm this law. The reactions we will use in this experiment are:

- (1) Solid sodium hydroxide dissolves in water to form an aqueous solution of ions.

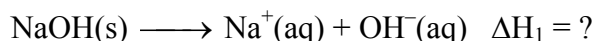

- (2) Solid sodium hydroxide reacts with aqueous hydrochloric acid to form water and an aqueous solution of sodium chloride.

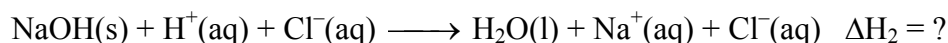

- (3) Solutions of aqueous sodium hydroxide and hydrochloric acid react to form water and aqueous sodium chloride.

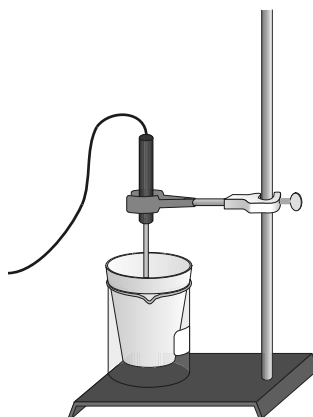
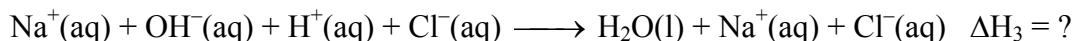


Figure 1

You will use a Styrofoam cup in a beaker as a calorimeter, as shown in Figure 1. For purposes of this experiment, you may assume that the heat loss to the calorimeter and the surrounding air is negligible. Even if heat is lost to either of these, it is a fairly constant factor in each part of the experiment, and has little effect on the final results.

OBJECTIVES

In this experiment, you will

- Combine equations for two reactions to obtain the equation for a third reaction.
- Use a calorimeter to measure the temperature change in each of three reactions.
- Calculate the heat of reaction, ΔH , for the three reactions.
- Use the results to confirm Hess's law.

MATERIALS

TI-Nspire handheld **or**
computer and TI-Nspire software
EasyTemp **or** Go!Temp, **or**
Temperature Probe and data-collection interface
50 mL of 1.0 M NaOH
50 mL of 1.0 M HCl
100 mL of 0.50 M HCl

100 mL of water
4.00 g of solid NaOH
ring stand
utility clamp
stirring rod
Styrofoam cup
250 mL beaker




PRE-LAB QUESTIONS

In the space below, combine two of the above equations algebraically to obtain the third equation. Indicate the number of each reaction on the shorter lines.

_____	_____
_____	_____
_____	_____

PROCEDURE

Reaction 1

1. Obtain and wear goggles.
2. Use a utility clamp to suspend a Temperature Probe from a ring stand as shown in Figure 1.
3. Connect the Temperature Probe to the data-collection interface. Connect the interface to the TI-Nspire handheld or computer. (If you are using an EasyTemp or Go!Temp, you do not need a data-collection interface.)
4. Choose New Experiment from the  Experiment menu. For this experiment, the default data-collection parameters for a temperature probe will be used (Rate: 2 samples per second; Duration: 180 seconds).
5. Place a Styrofoam cup into a 250 mL beaker as shown in Figure 1. Measure out 100.0 mL of water into the Styrofoam cup. Lower the Temperature Probe into the solution.
6. Weigh out about 2 g of solid sodium hydroxide, NaOH, and record the mass to the nearest 0.01 g in Table 2. Since sodium hydroxide readily picks up moisture from the air, it is necessary to weigh it and proceed to the next step without delay. **CAUTION:** Handle the NaOH and resulting solution with care. They are extremely corrosive.
7. Start data collection (). Monitor temperature (in °C) on the screen. It may take several seconds for the Temperature Probe to equilibrate at the temperature of the solution. After three or four readings at the same temperature have been obtained, add the solid NaOH to the Styrofoam cup. Using the stirring rod, stir continuously until the temperature has maximized and then begun to drop.
8. Data collection will stop after 3 minutes. If desired, you can stop data collection early (.

9. To confirm the initial (t_1) and maximum (t_2) temperature values, click any data point and use ► and ◀ move along the curve on the displayed graph. Record these values in Table 1.
10. Rinse and dry the Temperature Probe, Styrofoam cup, and stirring rod. Dispose of the solution as directed by your instructor.

Reaction 2

11. Click the Store Latest Data Set button (⌂) to save the data from Reaction 1. Repeat Steps 5–10, using 100.0 mL of 0.50 M hydrochloric acid, HCl, instead of water. Use the same amount of solid NaOH as before. **CAUTION:** *Handle the HCl solution and NaOH solid with care.*

Reaction 3

12. Click the Store Latest Data Set button (⌂) to save the data from Reaction 2. Repeat Steps 5–10, initially measuring out 50.0 mL of 1.0 M HCl (instead of water) into the Styrofoam calorimeter. In Step 5, instead of solid NaOH, measure 50.0 mL of 1.0 M NaOH solution into a graduated cylinder. After 3–4 temperature readings have been taken to determine the initial temperature of the 1.0 M HCl, add the 1.0 M NaOH solution to the Styrofoam cup. **CAUTION:** Handle the HCl and NaOH solutions with care.

DATA

Table 1			
	Reaction 1	Reaction 2	Reaction 3
Mass of solid NaOH	g	g	(no solid NaOH mass)
Final temperature, t_2	°C	°C	°C
Initial temperature, t_1	°C	°C	°C

PROCESSING THE DATA

1. Determine the mass, m , of 100 mL of solution for each reaction (assume the density of each solution is 1.00 g/mL).
2. Determine the temperature change, Δt , for each reaction.
3. Calculate the heat released by each reaction, q , by using the formula:

$$q = C_p \cdot m \cdot \Delta t \quad (C_p = 4.18 \text{ J/g}^\circ\text{C})$$

Convert joules to kJ in your final answer.

4. Find ΔH ($\Delta H = -q$).
5. Calculate moles of NaOH used in each reaction. In Reactions 1 and 2, this can be found from the mass of the NaOH. In Reaction 3, it can be found using the molarity, M , of the NaOH and its volume, in L.

