

Go!Link

Order Code: GO-LINK

Go! Link is a single-channel USB interface used to connect many Vernier sensors directly to a Windows® or Macintosh® computer. Some Vernier sensors can also be used when Go! Link is connected to a Chromebook™.




Note: Vernier products are designed for educational use. Our products are not designed, nor are they recommended, for any industrial, medical, or commercial process such as life support, patient diagnosis, control of a manufacturing process, or industrial testing of any kind.

Compatible Software

See www.vernier.com/product/golink for a list of software compatible with the Go! Link.

Getting Started

1. Connect Go! Link to a USB port and the sensor to the Go! Link.
2. Start the data-collection software. The sensor will be recognized and in most cases will display the live reading. **Note:** The thermometer display shown in the graph below is only displayed when using a temperature probe.
3. Click Collect to begin collecting data.

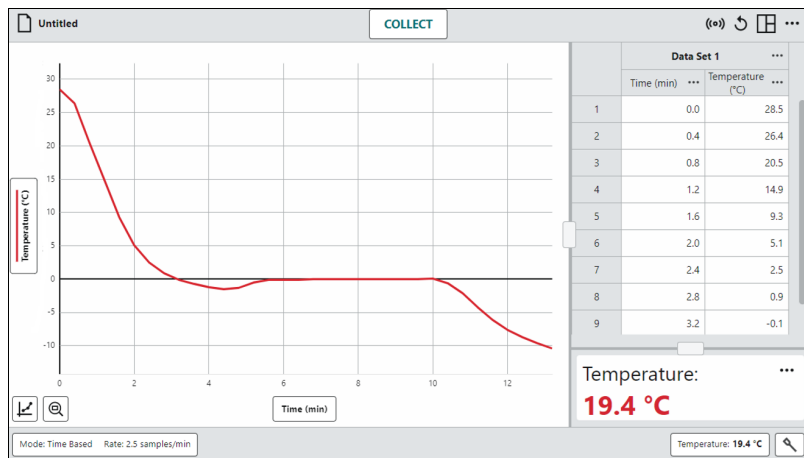

Specifications

USB specification	1.1
Maximum sample rate	200 samples per second
Resolution	12 bit

Accessories/ Replacements

Go! Link Teacher Pack (order code: GL-TP)

Save money by purchasing our Go! Link teacher pack. Use the 8 Go! Links to connect 8 sensors to 8 computers or connect multiple sensors to one computer or Chromebook.

Go! Temp® (order code: GO-TEMP)

Go!Temp is an easy-to-use temperature probe that connects directly to the USB port of a computer, LabQuest®, or Chromebook.

Go! Motion® (order code: GO-MOT)

Go! Motion is a motion detector that connects directly to the USB port of a computer, LabQuest, or Chromebook.

Troubleshooting

For troubleshooting tips and frequently asked questions regarding Go! Link, see www.vernier.com/til/2114

Warranty

Warranty information for this product can be found on the Support tab at www.vernier.com/go-link

General warranty information can be found at www.vernier.com/warranty



Vernier Science Education
13979 SW Millikan Way • Beaverton, OR 97005-2886
Toll Free (888) 837-6437 • (503) 277-2299 • Fax (503) 277-2440
info@vernier.com • www.vernier.com

Rev. 11/14/2024

Go! Link, Vernier LabQuest, Go! Motion, and other marks shown are our trademarks or registered trademarks in the United States.

All other marks not owned by us that appear herein are the property of their respective owners, who may or may not be affiliated with, connected to, or sponsored by us.

