

---

# Gas Pressure Sensor Bulb

Order Code GPS-BULB1, GPS-BULB4

## What's Included

- Gas Pressure Sensor Bulb (4 bulbs if GPS-BULB4 purchased)


## Using the Product

To measure grip strength, insert one of the stopper stems that comes with our gas pressure sensor into the exhaust end of the bulb. Then connect one of the tubing with a Luer-lock connector to the stopper stem and the other end to the gas pressure sensor.

## Tips

- Make sure the stopper stem is inserted securely and in the correct end of the bulb.
- Grasp the bulb with relaxed fingers. Do not curl fingers into the bulb.

**Note:** Vernier products are for educational use only.

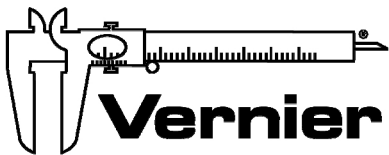

**Vernier Software & Technology**

13979 S. W. Millikan Way • Beaverton, OR 97005-2886

Toll Free (888) 837-6437 • (503) 277-2299 • Fax (503) 277-2440

info@vernier.com • www.vernier.com

Rev.12/16/19



Printed on recycled paper