

Exercise Heart Rate Strap

Order Code HR-STRAP



The Exercise Heart Rate Strap requires a Polar® Transmitter Module (HR-TRANS, not included). This combination is an excellent hands-free option for continuously monitoring heart rate before, during, and after exercise. Data are wirelessly transmitted to iPad®, LabQuest 2 or newer, or other mobile devices that are compatible with Bluetooth® wireless technology. For information on how to use the Exercise Heart Rate Strap with a Polar Transmitter Module, please see the Go Wireless® Exercise Heart Rate user manual at <http://www.vernier.com/gw-ehr>.

NOTE: Vernier products are designed for educational use. Our products are not designed nor are they recommended for any industrial, medical, or commercial process such as life support, patient diagnosis, control of a manufacturing process, or industrial testing of any kind.

Optional Accessories

Polar Transmitter Module (order code: HR-TRANS)

A module that transmits heart rate data wirelessly to Bluetooth Smart Ready devices or devices equipped to receive 5 kHz RF signals. For more information, see www.vernier.com/hr-trans.

Heart Rate Receiver (order code: HR-REC)

A 5 kHz RF receiver attached to a BTA connector. This allows the Polar Transmitter Module to communicate with Vernier interfaces that do not support Bluetooth Smart Ready devices. For more information, see www.vernier.com/hr-rec.

Disposal Instruction

When disposing of this electronic product, do not treat it as household waste. Its disposal is subject to regulations that vary by country and region. This item should be given to an applicable collection point for the recycling of electrical and electronic equipment. By ensuring that this product is disposed of correctly, you help prevent potential negative consequences on human health or on the environment. The recycling of materials will help to conserve natural resources. For more detailed information about recycling this product, contact your local city office or your disposal service.



The symbol, shown here, indicates that this product must not be disposed of in a standard waste container.

Exercise Heart Rate Strap

Order Code HR-STRAP



The Exercise Heart Rate Strap requires a Polar® Transmitter Module (HR-TRANS, not included). This combination is an excellent hands-free option for continuously monitoring heart rate before, during, and after exercise. Data are wirelessly transmitted to iPad®, LabQuest 2 or newer, or other mobile devices that are compatible with Bluetooth® wireless technology. For information on how to use the Exercise Heart Rate Strap with a Polar Transmitter Module, please see the Go Wireless® Exercise Heart Rate user manual at <http://www.vernier.com/gw-ehr>.

NOTE: Vernier products are designed for educational use. Our products are not designed nor are they recommended for any industrial, medical, or commercial process such as life support, patient diagnosis, control of a manufacturing process, or industrial testing of any kind.

Optional Accessories

Polar Transmitter Module (order code: HR-TRANS)

A module that transmits heart rate data wirelessly to Bluetooth Smart Ready devices or devices equipped to receive 5 kHz RF signals. For more information, see www.vernier.com/hr-trans.

Heart Rate Receiver (order code: HR-REC)

A 5 kHz RF receiver attached to a BTA connector. This allows the Polar Transmitter Module to communicate with Vernier interfaces that do not support Bluetooth Smart Ready devices. For more information, see www.vernier.com/hr-rec.

Disposal Instruction

When disposing of this electronic product, do not treat it as household waste. Its disposal is subject to regulations that vary by country and region. This item should be given to an applicable collection point for the recycling of electrical and electronic equipment. By ensuring that this product is disposed of correctly, you help prevent potential negative consequences on human health or on the environment. The recycling of materials will help to conserve natural resources. For more detailed information about recycling this product, contact your local city office or your disposal service.



The symbol, shown here, indicates that this product must not be disposed of in a standard waste container.

Warranty

Warranty information for this product can be found on the Support tab at www.vernier.com/gdx-acc

General warranty information can be found at www.vernier.com/warranty

Warranty

Warranty information for this product can be found on the Support tab at www.vernier.com/gdx-acc

General warranty information can be found at www.vernier.com/warranty



Vernier Science Education

13979 SW Millikan Way • Beaverton, OR 97005-2886
Toll Free (888) 837-6437 • (503) 277-2299 • Fax (503) 277-2440
info@vernier.com • www.vernier.com

Rev. 8/2/2024

Go Wireless is our trademark in the United States. All other marks not owned by us that appear herein are the property of their respective owners, who may or may not be affiliated with, connected to, or sponsored by us.



Vernier Science Education

13979 SW Millikan Way • Beaverton, OR 97005-2886
Toll Free (888) 837-6437 • (503) 277-2299 • Fax (503) 277-2440
info@vernier.com • www.vernier.com

Rev. 8/2/2024

Go Wireless is our trademark in the United States. All other marks not owned by us that appear herein are the property of their respective owners, who may or may not be affiliated with, connected to, or sponsored by us.

