

LabQuest[®] Charge Station

Order Code LQ3-CRG

The LabQuest Charge Station is used to charge LabQuest 3, LabQuest 2, LabQuest Stream[®], and original LabQuest interfaces.

For additional product information, visit www.vernier.com/lq3-crg

What's Included

- 4-slot Charge Station
- Charge Station Cable

Using the Product

Perform the following steps to charge your interface in the charge station:

1. Connect the power cable to the charge station and to an outlet.
2. Place the LabQuest interface in the slot so that its AC power port is easily accessible.
3. Remove the charge station connector cord from its storage hole and attach the connector to the AC port on the LabQuest interface.

Precautions and Operational Tips

- You can charge LabQuest interfaces for any amount of time, even for a few minutes between classes.
- The charge station is not water resistant and should not be exposed to a wet environment.
- LabQuest interfaces can be stored in the charger with the charge station connected to AC. LabQuest batteries cannot be overcharged.
- When not using a charging slot, place the connector cord in the storage hole. This protects the cord from damage.

Troubleshooting

The LabQuest interface charges with its individual AC adapter but not with the charge station.

- Make sure the charge station is connected to a working AC outlet.
- Attach the charge station connector to a LabQuest interface and verify that the unit is charging. For an original LabQuest, LabQuest 2, or LabQuest 3, you need to power the unit on and check the battery status icon.

Repair Information

If you have followed the troubleshooting steps and are still having trouble with your LabQuest Charge Station, contact Vernier Technical Support at support@vernier.com or call 888-837-6437. Support specialists will work with you to determine if the unit needs to be sent in for repair. At that time, a Return Merchandise Authorization (RMA) number will be issued and instructions will be communicated on how to return the unit for repair.

Accessories/Replacements

Replacement Parts

Charge Station Cable

Charge Station Cable Euro

Order Code

CRG-CB

CRG-CB-EURO

Disposal Instruction


When disposing of this electronic product, do not treat it as household waste. Its disposal is subject to regulations that vary by country and region. This item should be given to an applicable collection point for the recycling of electrical and electronic equipment. By ensuring that this product is disposed of correctly, you help prevent potential negative consequences on human health or on the environment. The recycling of materials will help to conserve natural resources. For more detail information about recycling this product, contact your local city office or your disposal service.



The symbol, shown here, indicates that this product must not be disposed of in a standard waste container.

Warranty

Vernier warrants this product to be free from defects in materials and workmanship for a period of five years from the date of shipment to the customer. This warranty does not cover damage to the product caused by abuse or improper use.

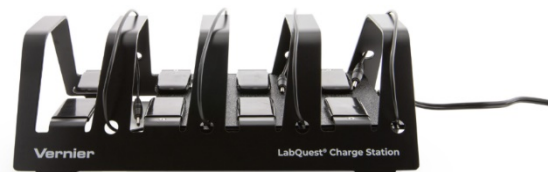



Vernier Software & Technology

13979 S. W. Millikan Way • Beaverton, OR 97005-2886

Toll Free (888) 837-6437 • (503) 277-2299 • FAX (503) 277-2440

info@vernier.com • www.vernier.com

Rev. 1/28/2021

Vernier LabQuest and other marks shown are our trademarks or registered trademarks in the United States.

All other marks not owned by us that appear herein are the property of their respective owners, who may or may not be affiliated with, connected to, or sponsored by us.