
Vernier Go Direct[®] Charge Station

Order Code GDX-CRG


PLEASE READ BEFORE USING PRODUCT!!

Please connect the included power supply to the GDX-CRG station and wall outlet **BEFORE** connecting any Vernier sensors.



Charging Instructions


Connect one or more supported Go Direct[®] devices or Go Wireless sensors to the Go Direct Charge Station. Charge the devices as recommended in each sensor's user manual.

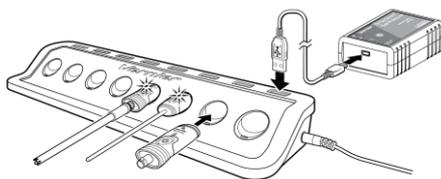

Charging Indicator

The charging indicator on each sensor is used to determine the charging status for that sensor. For charging LED behavior specific to your sensor, refer to the sensor's user manual.

Note: For wand-style sensors that fit into the oval holes in the charging station, verify each sensor shows its blue charging indicator light **when first connected** to the charge station (even if it is fully charged). If it does not, there may be an issue with the sensor. Contact Vernier's Support team for additional assistance.

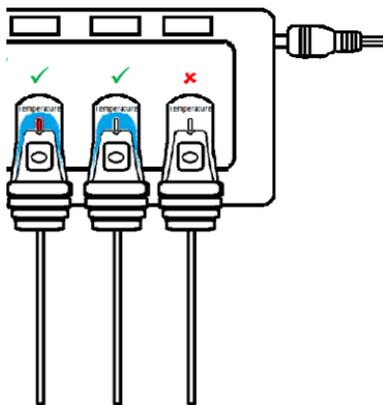

www.vernier.com/support/

1-888-837-6437

Precautions and Operational Tips

- You can charge devices for any amount of time, even for a few minutes between classes.
- Never insert metal objects into the charging slots.
- The Charge Station is not water resistant and should not be exposed to a wet environment.
- Do not insert a wet device into the Charge Station.
- Use only the power adapter provided with this unit.
- Do not attempt to charge a wand-style sensors in the oval charging slots if the sensor's blue charging-indicator light does not turn on at least briefly when first connected to the charging station that is connected to power. Doing so could shorten the life of your charging station.

Disposal Instruction

When disposing of this electronic product, do not treat it as household waste. Its disposal is subject to regulations that vary by country and region. This item should be given to an applicable collection point for the recycling of electrical and electronic equipment. By ensuring that this product is disposed of correctly, you help prevent potential negative consequences on human health or on the environment. The recycling of materials will help to conserve natural resources. For more detailed information about recycling this product, contact your local city office or your disposal service.



—The symbol, shown here, indicates that this product must not be disposed of in a standard waste container.

Warranty

Warranty information for this product can be found on the Support tab at www.vernier.com/gdx-crg

General warranty information can be found at www.vernier.com/warranty

Vernier Science Education
13979 SW Millikan Way • Beaverton, OR 97005-2886
Toll Free (888) 837-6437 • (503) 277-2299 • Fax (503) 277-2440
info@vernier.com • www.vernier.com

Rev. 2/12/2024

Go Direct and Go Wireless shown are our trademarks or registered trademarks in the United States. All other marks not owned by us that appear herein are the property of their respective owners, who may or may not be affiliated with, connected to, or sponsored by us.



Printed on recycled paper