

LabQuest Viewer Release Notes

Version 2.1

September 2026

System Requirements

This version of the LabQuest Viewer software is compatible with the following computer platforms:

- Macintosh — macOS 14, macOS 15, macOS 26
- Windows — Windows 10 (22H2 and newer), Windows 11

Users of older operating systems must continue using LabQuest Viewer v2.0 (macOS) and 2.0.1 (Windows).

Memory Requirements

LabQuest Viewer requires approximately 5.9 MB for macOS and 16 MB for Windows disk storage space.

Display Requirements

Computer/projector display resolution 1024x768 or higher is recommended. The minimum display resolution is 800x600.

Install Notes

LabQuest Viewer automatically discovers LabQuest devices on your local network using built-in operating system service discovery. No additional third-party networking software is required.

What's Changed in LabQuest Viewer 2.1

Macintosh

- Converted application to a Universal Binary to support newer macOS versions, including support for both Intel and Apple Silicon processors.
- Fixed issue where LabQuest screens were not displayed when added to My Lab.

Windows

- Converted to built-in Windows network service discovery (no separate client software required).

All Platforms

- Fixed issue where Check for Updates was causing an error.

LabQuest Viewer Change History

LabQuest Viewer 2.0.1 (*Windows only*)

- Fixed issue where LabQuest Viewer could leave LabQuest unresponsive to taps.

LabQuest Viewer 2.0

New Features

- New My Lab view displays live images of all LabQuest units in your lab.
- My Lab Setup allows you to add, arrange, and remove LabQuest devices in My Lab.
- Persistence of My Lab setup and software size/position values between sessions.
- Previous/Next buttons support quickly moving through LabQuest units while presenting.
- Select up to four LabQuest for side-by-side screen comparison.

Platform Support

- Macintosh — OS X 10.8, OS X 10.9, and OS X 10.10
- Windows — Windows 7 SP1, Windows 8, and Windows 8.1
- LabQuest Viewer makes use of Apple's Bonjour software for automatic device discovery.

LabQuest Viewer 1.1

Platform Support

- Windows — Windows XP, Windows Vista

LabQuest Viewer 1.0

Platform Support

- Macintosh — OS X 10.5 (Intel Macs only), OS X 10.6, and OS X 10.7